

Quadratech Instrument Cleaner

AGG373A

Application details

Ammonia free cleaner

A concentrated cleaning solution, suitable for the cleaning of instrument or watch parts. Formulated without ammonia to reduce fumes and odour in use. It is safe to all normal instrument materials and will remove tarnish, old oil and grease films, working quickly without causing damage.

Directions for use

Dilute one part of product with between 5 and 20 parts of water depending on the degree of soiling and tarnish present. Agitate the items to be cleaned in the solution for a few minutes, rinse thoroughly in clean water and dry. Best results will be obtained with ultrasonic agitation.

Please refer to product safety data sheet before use. Use only as directed.

Key product features

- NON AMMONIA FORMULATION
- ULTRASONIC COMPATIBLE
- SAFER IN USE